

# Outlet Cap & Gold Seal Assembly Tool

for Agilent 1100 HPLC Systems



**EASILY INSTALL  
GOLD SEAL INTO  
OUTLET CAP**



Put Outlet Cap on male part of Assembly Tool.



Push the Assembly Tool together.



Hold onto Outlet Cap and pull Assembly Tool apart.



Put Gold Seal over pin on male part of Assembly Tool.



Push the Assembly Tool together and press the Gold Seal in the Outlet Cap.



Pull the Assembly Tool apart and remove assembled Outlet Cap and Gold Seal.

**Description**

Outlet Cap and Gold Seal Assembly Tool for Agilent 1100 HPLC Systems

qty.

ea.

cat.#

24989

**Restek Replacement Parts for Agilent HPLC Systems**

• Meet or exceed OEM performance.

**Description**

Seal, Gold Disk (outlet)

Model #

1050 & 1100

Similar to Agilent part #

5001-3707

qty.

ea.

cat.#

25467

Outlet Cap

1050 & 1100

5062-2485

4-pk.

25139

Outlet Cap & Gold Seal Assembly

1050 & 1100

—

2-pk.

25140



**Chromatography Products**

Lit. Cat.# 580159

© 2007 Restek Corporation.

**HROM**alytic +61(0)3 9762 2034  
**ECH**nology Pty Ltd  
 Website NEW : www.chromalytic.com.au E-mail : info@chromtech.net.au Tel: 03 9762 2034 . . . in AUSTRALIA

**Australian Distributors**  
 Importers & Manufacturers  
 www.chromtech.net.au

