

IN THIS ISSUE

Editorial

Using Guard Columns and Retention Gaps in GC (Part 2) ..... 2

Environmental

Accurately Quantify PAHs Down to 5pg On-Column ..... 3

13 Minute Chlorophenoxyacid Herbicides Analysis ..... 6

Enhancing Air Monitoring Methods with Thermal Desorption ..... 8

Chemical/Petrochemical

Selecting a GC Column for Glycerin in Biodiesel ..... 10

Stable Sulfur & Mercury Sampling in Refineries ..... 12

Foods, Flavors & Fragrances

High Sensitivity Melamine |GC/MS Analysis of Cat Food ..... 14

Sample Preparation

Fast, Simple Sample Cleanup ..... 16

Pharmaceutical

Multi-task with an Ultra IBD Column ..... 18

Clinical/Forensics

Fast, Sensitive Analysis of Benzodiazepines by LC/MS/MS ..... 20

Accurate, Reproducible Amphetamines Analysis ..... 22

Tech Tip

Selecting the Right HPLC Guard Column ..... 26

HPLC Accessories

Waste Overflow Indicator for HPLC Systems ..... 27

GC Accessories

The Forgotten Septum ..... 24

Peak Performers FID Detector Replacement Parts ..... 28

Electron Multipliers for Mass Spectrometry ..... 30

Restek Trademarks

Allure, Alumaseal, Crossbond, Integra-Gap, Integra-Guard, MXT, Press-Tight, Rtx, Rxi, SeCure, Silcosteel, Siltek, Sulfinert, Uniliner, Restek logo.

Other Trademarks

Dacthal (Amvac Chemical Corp.), API 3200, Cliquid, Q Trap (Applied Biosystems/MDS SCIEX Instruments), Snap Seal (J.G. Finneran Associates, Inc.), SecureTD-Q (Markes International, UK), Parker (Parker Intangibles LCC Ltd.), Swagelok (Swagelok Company), Upchurch Scientific (Upchurch Scientific, Inc.), Valco (Valco Instruments Company, Inc.).

# Using Guard Columns and Retention Gaps in GC (Part 2)

Jaap de Zeeuw, International GC Consumables Specialist, Restek Corporation



Guard columns and retention gaps are used widely in gas chromatography (GC). Many users have difficulty understanding the difference between these two products, even though there is a significant difference in application. In Part 1 of this article we reviewed retention gaps, which mainly are used for focusing the sample components when introducing a large (liquid) sample directly onto the column. In contrast, guard columns are used to protect the analytical column from contamination. Guard columns and retention gaps both must be coupled to the analytical column, and this connection introduces a potential point of risk. A new approach is to integrate the retention gap directly into the analytical column tubing. By applying a “segment” coating technology the stationary phase can be deposited only in a certain part of the column allowing a deactivated section at the beginning. Column coupling is not required and maintenance is greatly simplified. Here we will review guard columns and discuss the new segment coating technology.

## Use of guard columns

The purpose of using guard columns is to protect the analytical column from contamination since the sample that is introduced is not always pure. Although the best chromatography is obtained with “clean” samples, the practical situation is that sample clean-up procedures are minimized and relative “dirty” samples are introduced onto the column. Samples can contain particulates, heavy components, derivatization reagents, ionic residues, acids, bases... all these compounds can interfere with the stationary phase and they will influence the separation process. Usually the degradation of column performance is a slow process but it will happen.

Most of the time the impurities accumulate in the first meter(s) of the column and by cutting off this section adequate separation is restored. Many users choose to connect a guard column in front of the analytical column. Such a guard column is deactivated and can be trimmed when contaminated and eventually replaced. Depending on the application, guard columns have a lifetime of 1 week up to 6 months. One has different choices for guard columns; a guard column can consist only of deactivated capillary, or it can be a coated capillary.

*Deactivated capillary tubing:* Deactivated fused silica tubing can be purchased by the meter and then a defined length can be coupled in front of the analytical column. Upon contamination, a section of the guard column is removed. When the whole guard is “consumed” a new guard column can be coupled. The disadvantage of cutting parts off of the guard column is that the column becomes shorter and this may affect retention times. However, if a similar length is always cut from the guard column, the change in retention time becomes very predictable. A deactivated guard column will also result in band focusing. If the injection is not optimal, there will be a focusing effect similar to that of a retention gap.

*Coated capillary tubing:* As the guard column needs to prevent contamination of the analytical column, a coated guard column can help as it has both the surface deactivation and also the stationary phase layer. The easiest and most economical way of using coated guard columns (or precolumns) is to buy two analytical columns. One we will use as a separation column and the second one will be used to make coated guard columns. From this second column we will cut 2m sections and couple a section in front of the analytical separation column. We can run our samples until contamination affects peak shape/response and then we can replace the guard with a new 2m section.

The system we have created will produce reproducible retention times as we always will replace the entire 2m coated guard column. Since the stationary phase is the same on the guard as on the analytical column, there will be no surprises. The coated guard column also will allow more aggressive samples/more contamination before it will give up. Lastly, we are able to cut 15 coated guard columns from a full 30m analytical column...that's also economical! However, if using a coated guard column, there will be no focusing effects.

Continued on page 31.