

Liner Cap Removing Tool for Thermo Scientific GCs:**Focus GC / TRACE™ GC / Ultra/TRACE™ GC x GC**

- Easily loosens the liner cap from the injector.
- Unique, ergonomic handle—easy to grip.



Remove septum cap, septum holder, septum, and septum support.



Place tool on liner cap. Align two pins on bottom of tool with two open slots on liner cap.



Turn counterclockwise to loosen liner cap.



Use tweezers (cat. #20101) to remove liner cap.

Description	Similar to TS part #	qty.	cat.#
Liner Cap Removing Tool for Thermo Scientific GCs	205 070 10	ea.	24937

Make Life Easier (MLE) Capillary Tool Kit for Thermo Scientific GCs**Includes:**

- capillary installation gauge for Thermo Scientific GCs
- liner cap removing tool for Thermo Scientific GCs
- 1/8", 3/16", and 1/4" nylon brushes
- 1/4", 3/8", and 3/16" stainless steel wire tube brushes
- stainless steel surface brush
- 6 stainless steel jet reamers (0.25–0.65mm OD)
- 1/4" x 5/16" open end wrench
- 3/8" x 7/16" open end wrench
- 6mm x 7mm open end wrench
- 8mm x 10mm open end wrench
- 16mm x 17mm open end wrench
- rubber-tipped slide-lock tweezers
- scoring wafers with handles
- inlet liner removal tool
- septum puller
- mini wool puller/insertor tool
- 4-inch tapered needle file
- swivel head flashlight
- mini hand drill set
- 15cm compact steel ruler
- pocket magnifier
- high temperature string (1 meter)
- pipe cleaner (12-inch)
- cotton tip swabs (pk. of 25)

Everything you need in one complete kit!



Description	qty.	cat.#
MLE Capillary Tool Kit for Thermo Scientific GCs	kit	22183

Transfer Line Reducing Kit for Thermo Scientific TRACE™ and Focus DSQ Mass Spectrometers

- Meets or exceeds manufacturer's performance.



Description	Similar to TS part #	qty.	cat.#
Transfer Line Reducing Kit for Thermo Scientific TRACE™ and Focus DSQ Mass Spectrometers	76458-2014s 76458-2009s A0101-03151	kit	22082