



No more burned fingers!

Inlet Liner Removal Tool

- Easily remove liner from injector—no more burned fingers.
- Made from high-temperature silicone.
- Won't chip or crack the liner.



Gently push the liner removal tool onto the liner in the injection port, with a slight circular motion.



Slowly pull the liner out of the injection port.



Use the liner removal tool to place a new liner into the injection port, avoiding hot metal surfaces.

Description

Inlet Liner Removal Tool

qty.

3-pk.

cat.#

20181

Inlet Liner Packing Tool

- Position wool reproducibly every time.
- Accurate to a specific, measured depth.

Eliminates user variation!



Loosen the nut on the side of the tool and adjust the gauge to the manufacturer's recommended depth.



Place a plug of loosely bound wool at the top of the inlet liner.



Insert the liner packing tool into the liner until the tool bottoms out. Remove the tool. The wool is now positioned correctly in the liner and the liner is ready for use.

Description

Inlet Liner Packing Tool

qty.


ea.

cat.#

20339

Mini Wool Puller/Inserter

Insert and remove wool plugs easily. Order a spare pack so you'll always have one available.

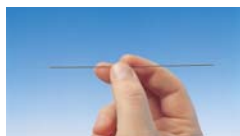

Description

Mini Wool Puller/Inserter

qty.

2-pk.

cat.#

20114

Rethreading Tool

- Repair worn or damaged threads.
- Multiple uses (fittings, injectors, detectors, etc.)
- Built-in guide to prevent cross-threading.

Achieve a better seal!



cat.# 23018



damaged threads

Repeated installation, removal, and temperature changes cause wear and damage to threads on GC parts. This can cause a poor seal, and oxygen can enter the system, compromising analytical results and possibly destroying expensive analytical columns.



Simply screw the rethreading tool onto the part, unscrew, and inspect the threads. Repeat as necessary. When done, wipe threads clean with methanol to remove any debris.



repaired threads

Description

Rethreading Tool for 1/16" compression fitting

qty.

ea.

cat.#

23016

Rethreading Tool for 1/8" compression fitting

ea.

23017

Rethreading Tool for 1/4" compression fittings (Agilent split/splitless injection ports)

ea.

23018

Rethreading Tool for 7/16" compression fitting (Varian injection ports)

ea.

23019

Rethreading Tool for 1/4" Varian-style capillary column fittings

ea.

21893