

## Outlet Cap and Gold Seal Assembly Tool for Agilent 1050 & 1100 HPLC Systems

Easily install the gold seal into the outlet cap.



Put Outlet Cap on male part of Assembly Tool.

Push the Assembly Tool together.

Hold onto Outlet Cap and pull Assembly Tool apart.

Put Gold Seal over pin on male part of Assembly Tool.

Push the Assembly Tool together and press the Gold Seal in the Outlet Cap.

Pull the Assembly Tool apart and remove assembled Outlet Cap and Gold Seal.



### Description

Description	qty.	cat.#
Outlet Cap and Gold Seal Assembly Tool for Agilent 1050 & 1100 HPLC Systems	ea.	24989
Seal, Gold Disk (outlet)	ea.	25467
Outlet Cap	4-pk.	25139
Outlet Cap & Gold Seal Assembly	2-pk.	25140



Outlet Cap & Gold Seal Assembly

## Restek Replacement Parts for Beckman HPLC Systems

Description	Model #	Similar to Beckman part #	qty.	cat.#
Inlet Check Valve Assembly	110A&B, 112, 114M, 116, 118, 125, 126, 127, 128 Pumps	240720	ea.	25439
Outlet Check Valve Assembly	110A&B, 112, 114M, 116, 118, 125, 126, 127, 128 Pumps	240721	ea.	25440
Inlet Check Valve Cartridge for Check Valve Assembly	110A&B, 112, 114M, 116, 118, 125, 126, 127, 128 Pumps	240620	ea.	25441
Outlet Check Valve Cartridge for Check Valve Assembly	110A&B, 112, 114M, 116, 118, 125, 126, 127, 128 Pumps	240621	ea.	25442
Guide Sleeve for Graphite Plunger Guide	110 Series Pumps	243713	ea.	25444
Outlet Check Filter Frit Assembly	110 Series Pumps	240619	ea.	25445
Piston Guide Assembly	110 Series and 112 Pumps	243045	ea.	25446
Plunger	110 Series and 112 Pumps	243053	ea.	25447
Plunger Seal	110 Series Pumps	887138	ea.	25448
Plunger Seal	112 Pumps	236797	ea.	25449
Sapphire Plunger	114M, 116, 118, 125, 126, 127, 128 Pumps	240714	ea.	25450
Plunger Seal	114M, 116, 118, 126, 127, 128 Pumps	241037	ea.	25451
Pump Seal, Gold*	114, 116, 125, 126, 127, 128 Pumps	241037	ea.	25452
Plunger Wash Seal	125, 126, 127, 128 Pumps	728772	ea.	25453
Deuterium Lamp	DU60, 62, 64, 65	596791	ea.	25454

\*Ultra-high molecular weight polyethylene (UHMWPE).



Outlet Check Valve Assembly



Sapphire Plunger

## Restek Replacement Parts for Hitachi HPLC Systems

Description	Model #	Similar to Hitachi part #	qty.	cat.#
SS Check Valve Cartridges (1/16" Ruby Ball)	655/6000/6200	AN0-085	2-pk.	25455
Inlet Check Valve Assembly	655/6000/6200	AN0-0833	ea.	25456
Outlet Check Valve Assembly	655/6000/6200	AN0-0834	ea.	25457
Plunger Assembly	655, 6000, 6200, 7100	810-1033	ea.	25458
Plunger Seal	L655A, 6000, 6200, 7100	655-1080	ea.	25459
Plunger Seal, Gold	655, 6000, 6200, 7100	655-1080	ea.	25460
Inlet Check Valve Assembly	L-7100	AN0-0836	ea.	25461
Outlet Check Valve Assembly	L-7100	AN0-0837	ea.	25462
Rotor Assembly for Dilutor Valve	AS-7200 Autosampler	810-3085	ea.	25463
Rotor Seal Kit	AS-7200, AS-7250 Injection Valve	AN0-0818	ea.	25464
Deuterium Lamp, Prealigned	L4000, L4200, L4250, L7400	885-3570	ea.	25465

