

restek
innovation!

Kit installs easily,
without special tools
or plumbing.




EZ No-Vent® GC Column-Mass Spectrometer Connector

for Varian Saturn 2000 Series Mass Spectrometers

- Change GC/MS columns in minutes without venting—100 µm transfer line throttles vacuum and prevents venting.
- Easy to install and maintain—no special tools or plumbing required.
- Gold-plated body for inertness.
- Deactivated transfer line keeps analytes focused; high-temperature polyimide ferrules eliminate leaks at the problematic transfer line fitting.
- Lower cost than other “no-vent” fittings.

We designed the EZ No-Vent® GC column-mass spectrometer connector to be simple and easy to use. A critical orifice in the EZ No-Vent® connector minimizes the amount of oxygen allowed into the MS source, eliminating the need for purge gas as is required for other manufacturer’s vent systems. This enables you to skip the lengthy vent and pump-down cycle otherwise required when you make a column change. This can save nearly a day of down-time with each column change. The EZ No-Vent® connector easily attaches to the MS source without special tools or extra plumbing.

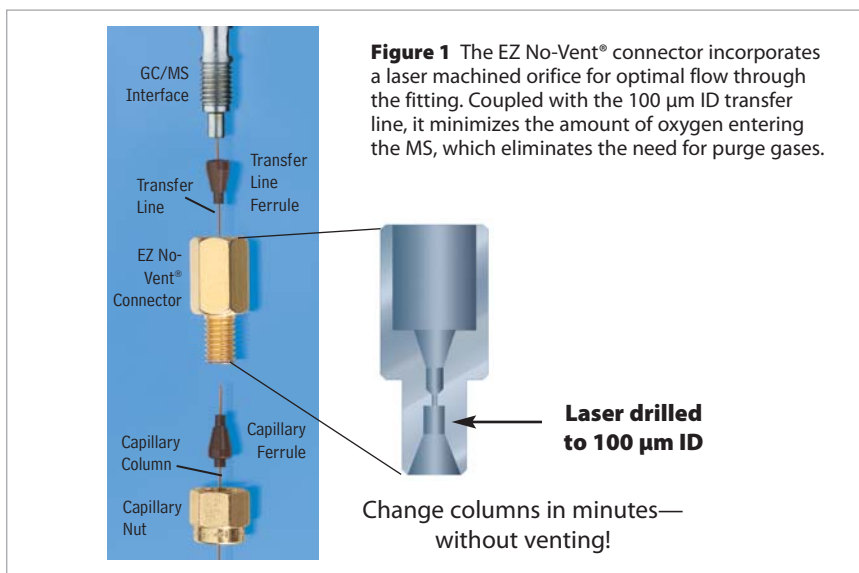

Description	qty.	cat.#	price
EZ No-Vent Connector Kit for Varian Saturn 2000 Series MSs			
Includes: EZ No-Vent Connector, two 0.4mm ID adaptor ferrules for capillary column, two 0.4mm ID ferrules for transfer line, 100µm deactivated transfer line (3 ft.), column plug, column nut	kit	22423	
Replacement ferrules for connecting capillary column to EZ No-Vent Connector:			
0.4mm ID (Polyimide)	2-pk.	21015	
0.5mm ID (Polyimide)	2-pk.	21016	
Replacement ferrules for connecting transfer line to EZ No-Vent Connector:			
0.4mm ID	2-pk.	21043	
Replacement 100µm deactivated transfer line	3 ft.	21018	
Replacement EZ No-Vent Column Nut	20-pk.	23100	
Replacement EZ No-Vent Plug	5-pk.	23112	
Open-End Wrenches, 1/4" x 5/16"	2-pk.	20110	