


Triglycerides

Triglycerides (Butter)

Rtx®-65TG

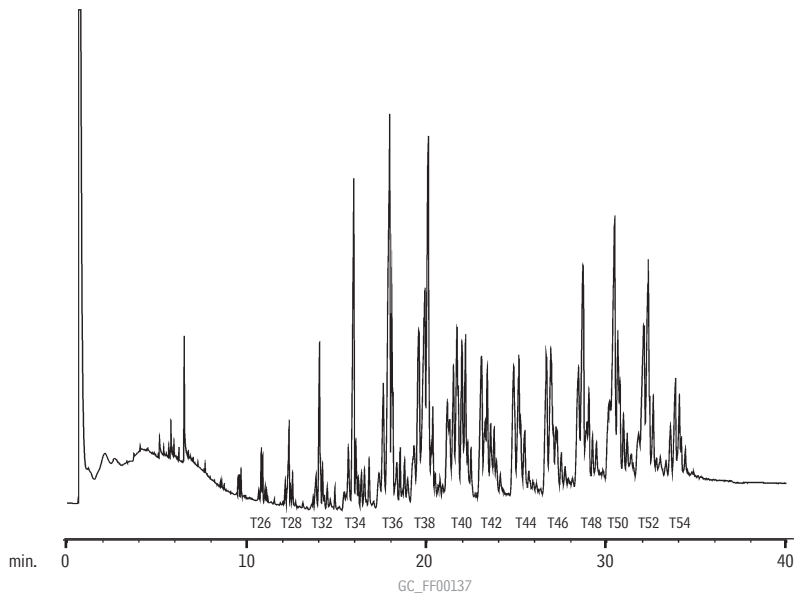

Column: Rtx®-65TG, 30m, 0.25mm ID, 0.10 μ m (cat.# 17008)
 Sample: 0.2 μ L cold on-column injection of 1% butterfat in isooctane
 Oven temp.: 80°C (hold 1 min.) to 240°C @ 30°C/min.
 to 360°C @ 4°C/min. (hold 5 min.)
 Det. temp.: 380°C
 Carrier gas: hydrogen
 Linear velocity: 70cm/sec.
 FID sensitivity: 16 x 10¹¹ AFS
 SPI injector: high performance capillary insert 60°C,
 300°C/min. to 400°C (hold 5 min.)



Al Carusone, Technical Service

Technical Service

Do you have a technical question? Restek's Technical Service group has answers! Drawing from our extensive libraries of technical information and many years of collective chromatography experience, the experts in Technical Service can help you from set-up to method development.

Contact us:

For quick answers to commonly asked questions any time of the day, visit www.restek.com/answers or contact us directly:

In the U.S.

Phone: 1-800-356-1688, ext. 4

Fax: 814-353-1568

e-mail: support@restek.com

Outside the U.S.

Contact your Restek representative.