

# Resprep™ Extraction Cells

for ASE® 200 Systems, Restek Enhanced Design

- Simpler design with fewer parts. Caps include frit, washer, PTFE O-ring, and threaded insert.
- Faster installation.
- Inner surfaces polished, for easier cleaning.
- Choose original equipment-equivalent stainless steel, or Siltek® deactivation for improved inertness.



**HROMalytic Chromatography**  
Australian Distributors **TECHnology** Products '08  
www.chromtech.net.au E-mail : info@chromtech.net.au Tel : +61 3 9762 2034 Fax : +61 3 9761 1169

**RESTEK**

**RESTEK**

**Chromatography Products**

www.restek.com 800-356-1688 • 814-353-1300

## Resprep™ Extraction Cells for ASE® 200 Systems, Restek Enhanced Design

These Restek replacement parts will keep your extraction system running smoothly. All parts are economically priced to save you money and are designed to meet or exceed the performance of the original manufacturer's parts.

To simplify replacing the frit and PEEK® washer in the ASE® 200 extraction cell cap, we have developed an enhanced cell design that eliminates the snap ring. The frit presses snugly into a groove in a new PEEK® washer design that fits into a recessed cavity in the enhanced extraction cell.



Description	Stainless Steel		Siltek® Treated	
	qty.	cat.#	qty.	cat.#
Extraction Cell Kit, Resprep™ for ASE® 200, 1mL	kit	25980	kit	25981
Extraction Cell Kit, Resprep™ for ASE® 200, 5mL	kit	25982	kit	25983
Extraction Cell Kit, Resprep™ for ASE® 200, 11mL	kit	25984	kit	25985
Extraction Cell Kit, Resprep™ for ASE® 200, 33mL	kit	25986	kit	25987
Extraction Cell Body, Resprep™ for ASE® 200, 1mL	ea.	25960	ea.	25961
Extraction Cell Body, Resprep™ for ASE® 200, 5mL	ea.	25962	ea.	25963
Extraction Cell Body, Resprep™ for ASE® 200, 11mL	ea.	25964	ea.	25965
Extraction Cell Body, Resprep™ for ASE® 200, 33mL	ea.	25966	ea.	25967
Extraction Cell Caps, Resprep™ for ASE® 200	2-pk.	25968	2-pk.	25969
PEEK® Seal/Frit Assembly, Resprep™ for ASE® 200	2-pk.	25970	2-pk.	25971
Frit, Resprep™ for ASE® 200	12-pk.	25972	12-pk.	25973

  

Description	qty.	cat.#
PEEK® Seal, Resprep™ for ASE® 200	12-pk.	25974
PEEK® Seal, Resprep™ for ASE® 200	48-pk.	25975
PTFE O-Rings for ASE® 200 & ASE® 300 Caps	100-pk.	26187
Viton® O-Rings for ASE® 200 & ASE® 300 Caps	50-pk.	26188



## 2-in-1 Filter/O-Ring Insertion Tool Kit for ASE® 100/200/300

Inserting a Filter (ASE® 100/200/300 systems), using Filter Insertion Attachments on Resprep™ Tool Handle.



Screw the appropriate size attachment onto the end of the Resprep™ Tool Handle.



Place a filter at the top of the extraction cell.



Push the filter to the bottom of the extraction cell.



2-in-1 Filter/O-Ring Insertion Tool Kit

Inserting an O-Ring (ASE® 100/200/300 systems), using Resprep™ Tool Handle.



Place the O-ring over the tip of the tool.



Insert the tool into the center hole of the extraction cell cap.



Press the tool firmly inside the cap until the O-ring snaps into place.



Resprep™ Tool Handle

Description	qty.	cat.#
2-in-1 Filter/O-Ring Insertion Tool Kit for ASE® 100/200/300 (includes Resprep™ Tool Handle and Filter Insertion Attachments)	kit	26181
Resprep™ Tool Handle for ASE® 100/200/300	ea.	26182
Filter Insertion Attachments for ASE® 100/200/300 (1mL, 5mL, 11mL, 33mL)	4-piece set	26183



Filter Insertion Attachments

Restek Trademarks:  
Resprep, Siltek, Restek logo.

Other Trademarks:  
ASE (Dionex Corporation), PEEK (Victrex plc), Viton (E.I. du Pont de Nemours & Co., Inc.)



Lit. Cat.# 580140

© 2007 Restek Corporation.



www.chromtech.net.au E-mail : info@chromtech.net.au Tel : +61 3 9762 2034 Fax : +61 3 9761 1169

