

Minnesota Department of Agriculture List 1 Pesticides

Minnesota Ag List 1 Pesticides Mix A

(16 components)

acetochlor	metolachlor
alachlor	metribuzin
atrazine	pendimethalin
cyazazine	prometon
desethylatrazine	propachlor
desisopropylatrazine	propazine
dimethenamid*	simazine
ethalfluralin	trifluralin

200ppm each in acetone, 1mL/ampul

cat. # 32406 (ea.)

Minnesota Ag List 1 Pesticides Mix B

chlorpyrifos	phorate
EPTC	terbufos
fonofos	triallate

200ppm each in acetone, 1mL/ampul

cat. # 32407 (ea.)



Jonathan Keim

Technical Service

Supervisor


7+ years of service!

*Added to Minnesota Department of Agriculture List 1 pesticide (neutrals) incident investigation requirements, effective January 1, 2000.¹ CAS # 87674-68-8 manufactured by several companies under various trade names.

¹Analytical Lists for Pesticide Incident Investigations, Minnesota Department of Agriculture, Guidance Document 26 (3/99), St. Paul, MN. For a copy, visit their web site at: www.mda.state.mn.us

Minnesota Dept. of Agriculture List 1 Pesticides

- | | | | |
|---------------------------------|----------------------|-------------------|--|
| 1. EPTC**
(eptam, eradicane) | 6. desethylatrazine* | 12. terbufos** | 18. alachlor* |
| 2. propachlor* | 7. phorate** | 13. fonofos** | 19. metolachlor* |
| 3. ethalfluralin* | 8. prometon* | 14. triallate** | 20. chlorpyrifos**
(chlorpyrifos ethyl) |
| 4. trifluralin* | 9. simazine* | 15. dimethenamid* | 21. cyanazine* |
| 5. desisopropylatrazine* | 10. atrazine* | 16. acetochlor* | 22. pendimethalin* |
| | 11. propazine* | 17. metribuzin* | |

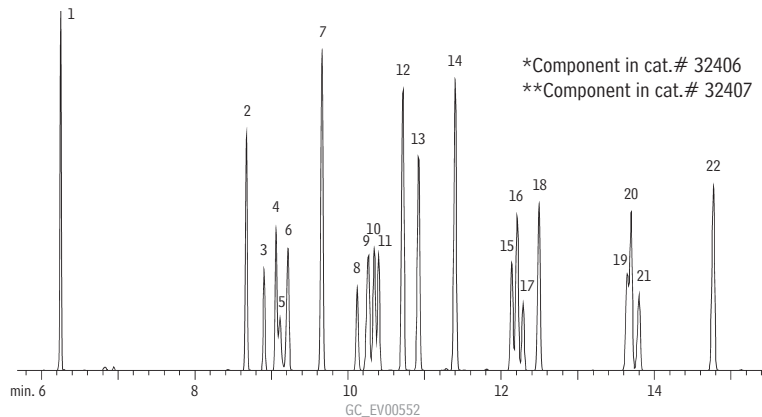

Column: Rtx®-5SiIMS, 30m, 0.28mmID, 0.5µm (cat.# 12794)
 Inj.: 1µL splitless 0.5 min., 10ppm calibration standard in methylene chloride/acetone, 10ng on column concentration
 Oven temp.: 35°C (hold 0.5 min.) to 160°C @ 50°C/min. to 190°C @ 10°C/min. to 300°C @ 5°C/min.
 Inj. temp: 250°C
 Det: Agilent 5973 MS, scan 35-550amu

free data

Available on Our Website: Lot Certificates, Data Packs, and MSDSs

For complete information detailing manufacturing and testing for Restek inventoried reference standards, visit our website at www.restek.com.

To view lot certificates and/or an MSDS, enter the catalog number of the product in the Search feature. For a free data pack (Adobe® PDF file), enter the catalog number and lot number of the product.