



Raspberry Flavor - Rt™-βDEXsa

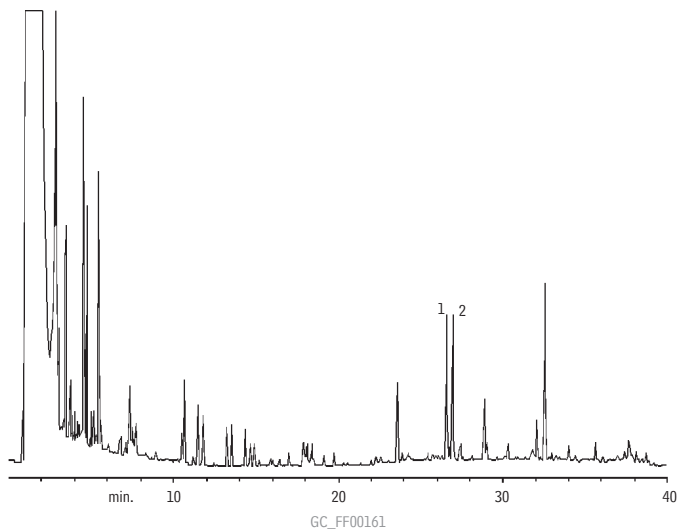

1. (R)-α-ionone
2. (S)-α-ionone

Column: Rt™-βDEXsa, 30m, 0.32mm ID, 0.25μm (cat.# 13108)
 Oven temp.: 60°C (hold 2 min.) to 200°C @ 3°C/min.
 Carrier gas: helium, 60cm/sec. set @ 60°C
 Det.: MS @ 220°C


Peach/Vanilla Flavor - Rt™-βDEXsa

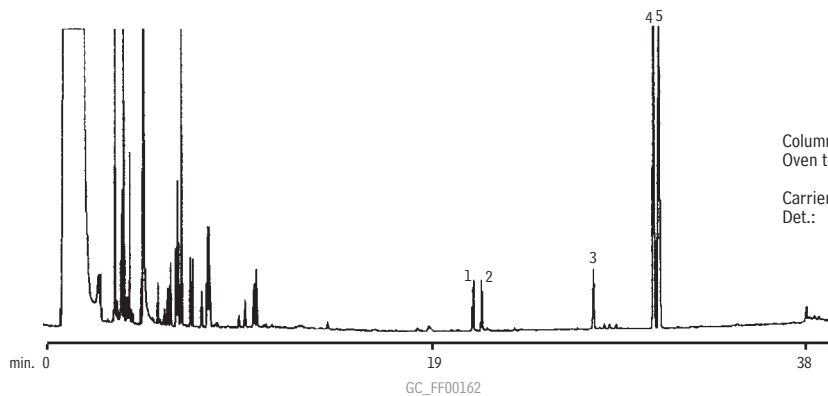

1. (R)-γ-octalactone
2. (S)-γ-octalactone
3. (R)-γ-decalactone
4. (R)-γ-undecalactone
5. (S)-γ-undecalactone

Column: Rt™-βDEXsa, 30m, 0.32mm ID, 0.25μm (cat.# 13108)
 Oven temp.: 60°C (hold 2 min.) to 100°C @ 15°C/min.
 to 220°C @ 3°C/min.
 Carrier gas: helium, 60cm/sec. set @ 60°C
 Det.: MS @ 220°C

Bergamot Flavor - Rt™-βDEXse

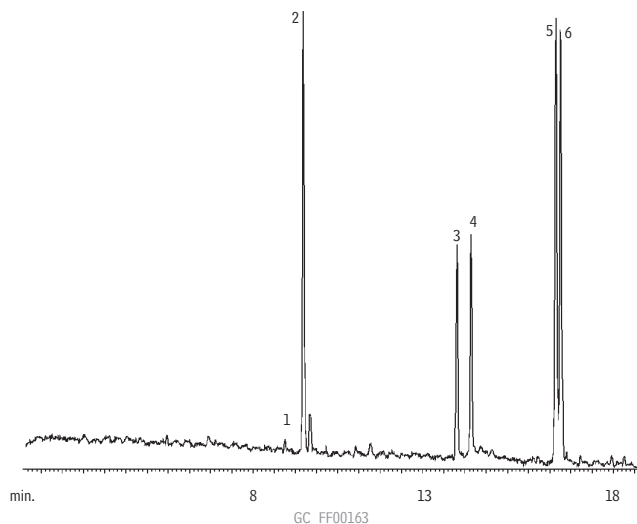

1. (S)-limonene
2. (R)-limonene
3. (R)-linalool
4. (S)-linalool
5. (R)-linalyl acetate
6. (S)-linalyl acetate

Column: Rt™-βDEXse, 30m, 0.32mm ID, 0.25μm (cat.# 13106)
 Oven temp.: 40°C (hold 1 min.) to 200°C @ 4°C/min.
 Carrier gas: helium, 60cm/sec. set @ 40°C
 Det.: MS @ 220°C