



Abused Inhalants

Rtx®-BAC1 & Rtx®-BAC2

Rtx®-BAC1

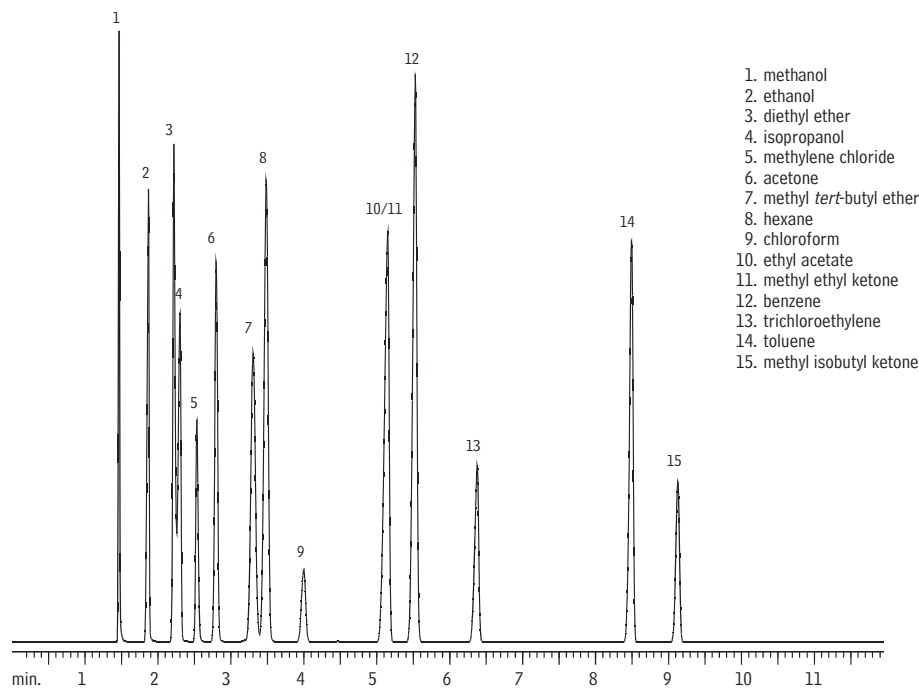

Rtx®-BAC2

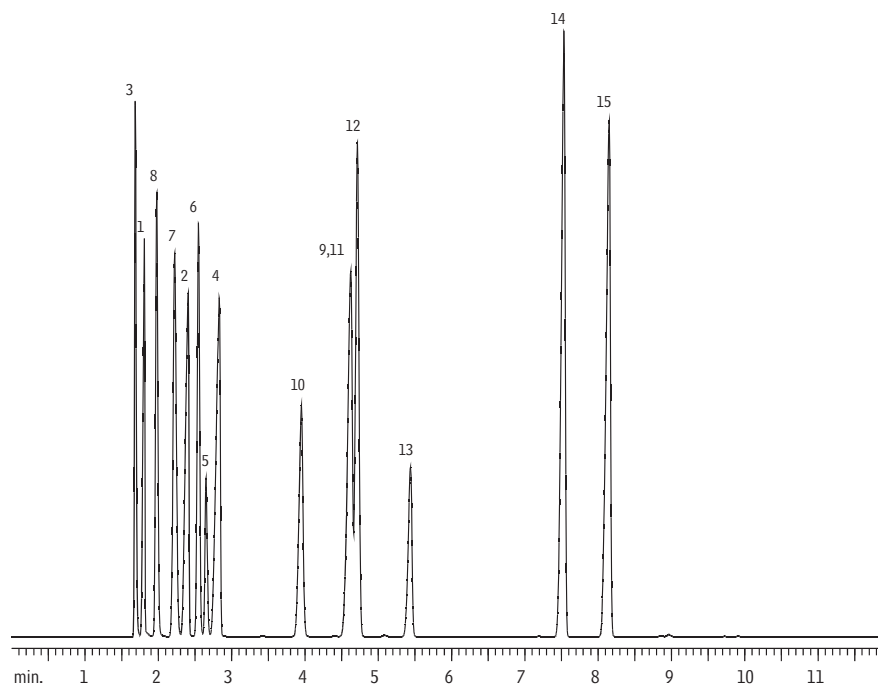

GC_PH00446

Columns: Rtx®-BAC1, 30m, 0.32mm ID, 1.80 μ m (cat.# 18003)
Rtx®-BAC2, 30m, 0.32mm ID, 1.20 μ m (cat.# 18002)

Oven temp.: 40°C (hold 4 min.) to 120°C @ 10°C/min.

Carrier gas: helium

Linear velocity: 50cm/sec.

Det.: FID

Sample: 250 μ L headspace

Transfer line temp.: 125°C

Sample loop temp.: 125°C

Vial temp.: 70°C

Acknowledgement: Analyses performed using a Tekmar 7000 Headspace Autosampler on loan courtesy of Tekmar-Dohrmann.