

Aromatics Analysis

D3606 Application Column (2 column set)

- Complete separation of ethanol and benzene, with a resolution value > 3.00.
- Accurate quantification of benzene and toluene.
- Fully conditioned two column set—ready to use out of the box.
- A chromatogram is provided with each column set demonstrating conformance to the revised ASTM method.

free literature

Resolve Benzene and Toluene in Spark Ignition Fuels Containing Ethanol

Download your free copy from

www.restek.com

lit. cat.# 580227




Conforms to the specifications established in ASTM method D3606-07 for the quantitation of benzene and toluene in spark ignition fuel containing ethanol.

Description	cat.#*	price
D3606 Application Column (2 column set)**		
Column 1: 6' (1.8m), 1/8" OD, 2.0mm ID, nonpolar Rtx-1		
Column 2: 16' (4.9m), 1/8" OD, 2.0mm ID, proprietary packing material	83606-	

*Please add column instrument configuration suffix number to cat.# when ordering. See page 143.

**The column set is designed to accommodate both valve injection and/or syringe injection. Column 1 is configured with a 2" inlet void to facilitate on-column injection. The inlet is identified on both column 1 and column 2. Note: The inlet of column 2 is identified for proper orientation for connection to the valve.

Gasoline containing ethanol on a D3606 Application Column set.

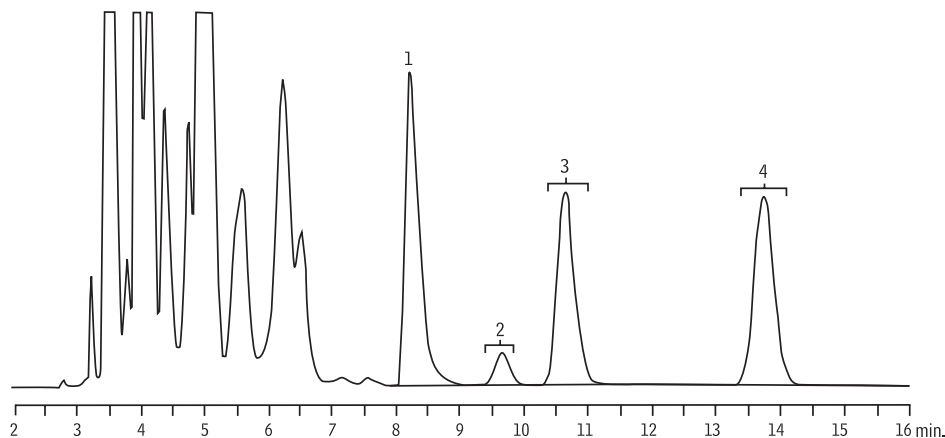

1. ethanol
2. benzene
3. 2-butanol
4. toluene

GC_PC01079

Column: D3606 Application Column (2 column set, cat.# 83606-800)
 Column 1: nonpolar Rtx®-1, 6' (1.8m), 1/8" OD, 2.0mm ID
 Column 2: proprietary packing material, 16' (4.9m), 1/8" OD, 2.0mm ID
 Sample: 1.5µL gasoline with internal standard
 Inj.: 200°C
 Backflush: 3 min.
 Carrier gas: helium, constant flow
 Flow rate: 20mL/min.
 Oven temp.: 135°C, isothermal
 Det.: TCD @ 200°C

Chromatogram courtesy of Boguslaw Dudek, Conoco Phillips, Linden, NJ.