

restek innovation!

No more burned fingers!

Inlet Liner Removal Tool

- Easily remove liner from injector—no more burned fingers.
- Made from high-temperature silicone.
- Won't chip or crack the liner.



20181



Gently push the liner removal tool onto the liner in the injection port, with a slight circular motion.



Slowly pull the liner out of the injection port.



Use the liner removal tool to place a new liner into the injection port, avoiding hot metal surfaces.

Description	qty.	cat.#	price
Inlet Liner Removal Tool	3-pk.	20181	




20339

Inlet Liner Packing Tool

- Position wool reproducibly every time.
- Accurate to a specific, measured depth.
- Can be used with all manufacturer's liners.



Loosen the nut on the side of the tool and adjust the gauge to the manufacturer's recommended depth.

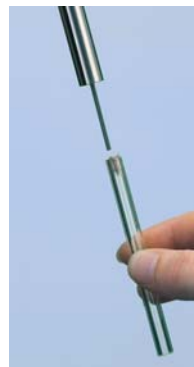

Place a plug of loosely bound wool at the top of the inlet liner.



Insert the liner packing tool into the liner until the tool bottoms out. Remove the tool. The wool is now positioned correctly in the liner and the liner is ready for use.

Eliminates user variation!

Description	qty.	cat.#	price
Inlet Liner Packing Tool	ea.	20339	

No more burned fingers!



26262

26261


The Claw and The Claw Holder Kit

- Easily removes hot liners from injection ports.
- 4mL vials (not included) can be replaced when dirty.

Never again will you have to burn your fingers removing a hot injection port liner. The Claw safely and cleanly removes liners, O-rings, or other small objects from the injection port. You can then place the hot objects in a clean 4 mL vial situated in the Claw holder until ready for reuse.



Description	qty.	cat.#	price
The Claw	ea.	26261	
The Claw Holder Kit (includes The Claw and holder)	kit	26262	
WISP 48 Snap Seal Vial	100-pk.	24658	

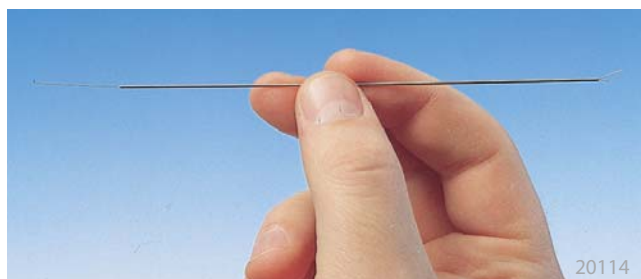

20114

Mini Wool Puller/Inserter

Insert and remove wool plugs easily. Order a spare pack so you'll always have one available.

Description	qty.	cat.#	price
Mini Wool Puller/Inserter	2-pk.	20114	

Sky

Introducing Sky™ Liners

Inlet Liners

See page 206 for more details.