

**Septum Nut Removal Tool**

for Agilent 5890/6890/6850/7890 GCs

- Easily remove the septum nut without touching the heated nut—no more burned fingers!
- Unique, ergonomic handle—easy to grip.
- Nut remains in tool for quick reattachment.

restek  
**innovation!**



24918



Slip tool over septum nut...



loosen nut...



and remove, avoiding hot metal surfaces.



Septum nut remains in tool until reinstalled.

**Liner Cap Removing Tool**

for Thermo Scientific GCs: Focus GC / TRACE GC / Ultra/TRACE GC x GC

- Easily loosens the liner cap from the injector.
- Unique, ergonomic handle—easy to grip.

restek  
**innovation!**



24937



Remove septum cap, septum holder, septum, and septum support.



Place tool on liner cap. Align two pins on bottom of tool with two open slots on liner cap.



Turn counter-clockwise to loosen liner cap.



Use tweezers (cat. #20101) to remove liner cap.

Description	qty.	cat.#	price
Septum Nut Removal Tool for Agilent 5890/6890/6850/7890 GCs	ea.	24918	

Description	Similar to TS part #	qty.	cat.#	price
Liner Cap Removing Tool for Thermo Scientific GCs	205 070 10	ea.	24937	





**Complete Product Offering**  
to support AOAC (2007.01) & European (EN 15662) QuEChERS Methods

- Comprehensive portfolio of QuEChERS tubes.
- Method-specific internal and QC standards.
- Rxi®-5Sil MS Capillary column
- Ultra Aqueous C18 LC Column.
- Q-sep™ 3000 Centrifuge

[www.restek.com/quechers](http://www.restek.com/quechers)