




25325

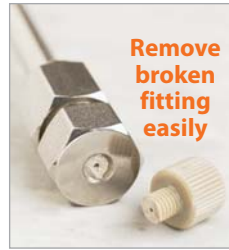

Remove broken fitting easily

Use the flat side of the Piston Seal Insertion Tool to seat a Waters face seal.



PEEK™ Fitting Extractor

Drill into the broken fitting, then screw the extractor into the fitting and remove it easily. Eliminates the need for heat or other techniques that could damage your column.



Description	qty.	cat.#	price
PEEK Fitting Extractor	ea.	25325	

HPLC Piston Seal Insertion Tool

Simplify pump maintenance: use one end to remove your old seal, then simply slip your new seal on the other end and push it flush into position. The tool cannot mar the surrounding metal surface of the pump housing.



Remove a seal



Seat a seal



25321

ValvTool Wrench

The ValvTool is a time-saving device that provides easy access to many hard-to-reach Rheodyne® or Valco® valves. For 1/4-inch nuts.

Description	qty.	cat.#	price
ValvTool Wrench	ea.	25321	

Description	qty.	cat.#	price
HPLC Piston Seal Insertion Tool	ea.	21356	

Outlet Cap and Gold Seal Assembly Tool for Agilent 1050 & 1100 HPLC Systems

Easily install the gold seal into the outlet cap.



Put Outlet Cap on male part of Assembly Tool.

Push the Assembly Tool together, then hold on to Outlet Cap and pull Assembly Tool apart.

Put Gold Seal over pin on male part of Assembly Tool.

Push the Assembly Tool together and press the Gold Seal into the Outlet Cap.

Pull the Assembly Tool apart and remove assembled Outlet Cap and Gold Seal.



25140

Description	qty.	cat.#	price
Outlet Cap and Gold Seal Assembly Tool for Agilent 1050 & 1100 HPLC Systems	ea.	24989	
Seal, Gold Disk (outlet)	ea.	25467	
Outlet Cap	4-pk.	25139	
Outlet Cap & Gold Seal Assembly	2-pk.	25140	