

**Resprep® Tools** for ASE® Systems

- Use to insert filter in extraction cell or O-ring in cell cap.
- Fits all extraction cells, except 1 mL size.

**2-in-1 Filter/O-Ring Insertion Tool Kit for ASE® 100/200/300**

Inserting a Filter (ASE® 100/200/300 systems)

Using Filter Insertion Attachments on Resprep® Tool Handle.




Screw the appropriate size attachment onto the end of the Resprep® Tool Handle.



Place a filter at the top of the extraction cell.



Push the filter to the bottom of the extraction cell.

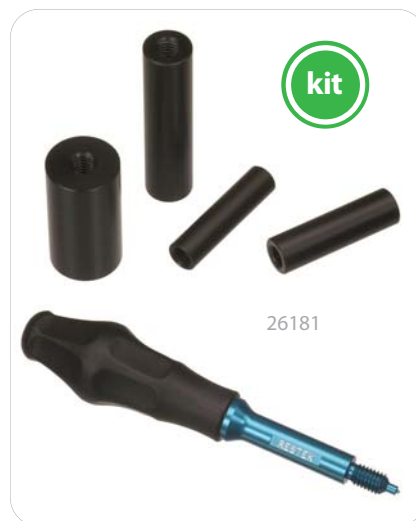

26181

Inserting an O-Ring (ASE® 100/200/300 systems)

Using Resprep® Tool Handle.



Place the O-ring over the tip of the tool.



Insert the tool into the center hole of the extraction cell cap.



Press the tool firmly inside the cap until the O-ring snaps into place.



26182



26183

Description	qty.	cat.#	price
2-in-1 Filter/O-Ring Insertion Tool Kit for ASE 100/200/300 (includes Resprep Tool Handle and Filter Insertion Attachments)	kit	26181	
Resprep Tool Handle for ASE 100/200/300	ea.	26182	
Filter Insertion Attachments (1mL, 5mL, 11mL, 33mL) for ASE 100/200/300	4-piece set	26183	

**Retaining Ring Pliers** for ASE® 100/200/300 Systems

- Can be used for internal or external retaining rings.
- Works to remove retaining ring in all ASE® cell caps.

Description	qty.	cat.#	price
Retaining Ring Pliers for ASE 100/200/300	ea.	26185	



26185

**Cell Cleaning Brushes** for ASE® 100/200/300 Cells

- Firm bristle brushes for easy cell cleaning and removal of solid samples.
- Range of sizes to fit all extraction cells for ASE® 100, 200, and 300.

Description	qty.	cat.#	price
Cell Cleaning Brushes	3-pk.	23999	

