

## Air Actuators

Air actuators offer reliable performance under the most stringent conditions. Low gas consumption and lightweight, compact construction make the air actuator suitable for aerospace flight hardware applications as well as laboratory or process applications.

The standard air actuator is rated for up to 80 psig at temperatures up to 70°C. Generally speaking, valves which will be heated require a standoff assembly, which locates the air actuator out of the heated zone and supports both the valve and actuator. A high temperature model permits both valve and actuator to be mounted within an oven (175°C maximum), but it is not recommended for use below 50°C.

### Air Actuators for Selectors

The recommended method for implementing a selector (multi-position) air actuator requires only a single 4-way solenoid. Up to 80 psig may be used without damaging the valve or actuator. Bottled instrument air or nitrogen is recommended.

If plant air from compressors must be used, an oil separator and water dryer are required.

Multiposition air actuators include a rotary switch which may be connected to a digital readout or your own design.




### Standard air actuators

*for selectors*

Temperature range 0-70°C  
Standoff version includes a 2" standoff. 3", 4", and 6" standoffs are also available.

Description	With closemount assembly	With 2" standoff assembly	With standoff mounting hardware
	Prod No	Prod No	Prod No
4 position	A4	A42	A4S
6 position	A6	A62	A6S
8 position	A8	A82	A8S
10 position	A10	A102	A10S
12 position	A12	A122	A12S
16 position	A16	A162	A16S

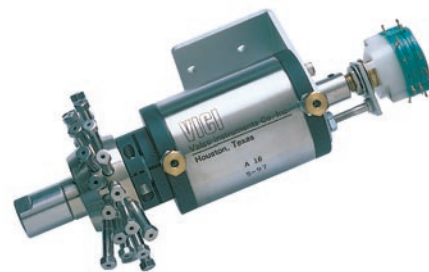

### High temperature air actuators

*for selectors*

Temperature range 50-175°C  
Standoff version includes a 4" standoff. 2", 3", and 6" standoffs are also available.

Description	With closemount assembly	With 4" standoff assembly	With standoff mounting hardware
	Prod No	Prod No	Prod No
4 position	AT4	AT44	AT4S
6 position	AT6	AT64	AT6S
8 position	AT8	AT84	AT8S
10 position	AT10	AT104	AT10S
12 position	AT12	AT124	AT12S
16 position	AT16	AT164	AT16S

#### TECH TIP

The actuator's rotation must be properly matched to the valve's. If you are converting a manual valve to air actuation and have any doubts about which actuator and hardware you need, call our sales or technical staff for assistance.

#### MORE INFORMATION

PFAF . . . . . page 199  
Position feedback

#### Mounting Hardware

Closemount hardware . . . . . page 208  
Right angle drive . . . . . 204  
Standoff assembly . . . . . 205  
Standoff mounting hardware . . . . . 205

#### ORDER TIP

To purchase a **valve with an air actuator installed**, go directly to valve ordering information.

### Replacement O-rings

Includes a complete set of O-rings for a multiposition air actuator.

Description	Prod No	Price
Standard	ORMP	\$16
High temp	ORTMP	20

