

Bench Top XL Crimper : *Crimper heads NOT Included > Order Separately !*

WARNING *see PDF www.chromtech.net.au/pdf2/XL BenchCrimper -Warning.pdf*

Take CARE in adjusting height of the Crimp Head "Holder" as it is relatively heavy (a few kg) AND needs Hand Support !

Otherwise it can inadvertently "crash" to the base and cause damage !

TO . . .

the Crimp Head jaws themselves

OR

to the base plate !

- *the "black-painted" Mild Steel base may be become scratched.*
 - *particularly if the head jaws are allowed to touch the base plate and the head rotated at the same time > easily causing an unsightly scratch.*
 - *normally the painted surface is fairly hardy but should be protected from use with aggressive solvents.*
- *a Granite Base can even crack !*
- *Both the SS and Granite base versions are also easy to clean > some solvent resistance eg alcohol / IPA wipe etc.*
- *Use the Stainless steel base version if autoclaving and sterility is required.*

ALWAYS use the "foam" protector between the plate and crimper jaws when the device is in transit or when the Bench Crimper is NOT being used !

> a strip of cardboard is a suitable alternative

Maintenance ?

None normally required ! apart from cap height/pressure adjustment

Except (maybe) in a Sterile dust-free room > very very light oil lubrication of jaws might be required if jaws tend to stick after extensive use !

