

Prevent Dangerous Spills with the Jour Level-Minder

- ◆ **Monitors Waste-Bottle Level - Audible and Visible Alarm Indicates 'Full' Condition Before Spills Occur**
- ◆ **Attaches to the Top of Standard Solvent Bottles with CL-45 threads**
- ◆ **Equipped with Two Convenient 'Waste' Connections**
- ◆ **Safe, Low-Voltage, Battery Operates Power Supply**

Anyone who has ever cleaned-up an overflowing HPLC waste bottle will appreciate Jour's Level-Minder™, the convenient waste-bottle level monitor that tells you, audibly and visibly, when your waste bottle is nearing full condition. Messy and dangerous spills are prevented, avoiding unnecessary hazardous waste exposure.

Level-Minder™ sits on top of the waste bottle with an inert float inside the bottle's neck. The self-contained, inherently safe, battery-operated system senses the float's position. When the float is at its lowest level, indicating the absence of solvent near the top of the bottle, Level-Minder™ green LED flashes, letting you know **Level-Minder™** is functioning and that the waste-bottle is OK. As the waste solvent reaches the float, **Level-Minder™** senses

the float's movement, activating a flashing red LED and audible alarm. The alarm activates well before the bottle overflows, giving you ample time to empty or change the bottle.



Level-Minder™ is self-contained requiring no additional bench space, wiring or fluid lines. Level-Minder™ fits on any bottle with CL-45 threads, the standard for most solvent bottles. Purchase the **Level-Minder™** alone, or with a safety-coated solvent bottle. Two convenient 1/4-28 TPI female ports accept waste solvent via 1/1 6" or 1/8" tubing. Flange-Free™ fittings for these connections are included. Two 'AA' batteries (included) power **Level-Minder™** for months of routine operation.

Why risk the danger, inconvenience, and mess from waste-bottle overflow? Add **Level-Minder™** to all your LC systems.

A. When the float is at its lowest level, the green LED flashes.

B. As solvent reaches the float the alarm is activated and the red LED flashes.

C. Two 1/4-28 TPI ports accept waste solvents via 1/1 6" or 1/1 8" OD tubing.

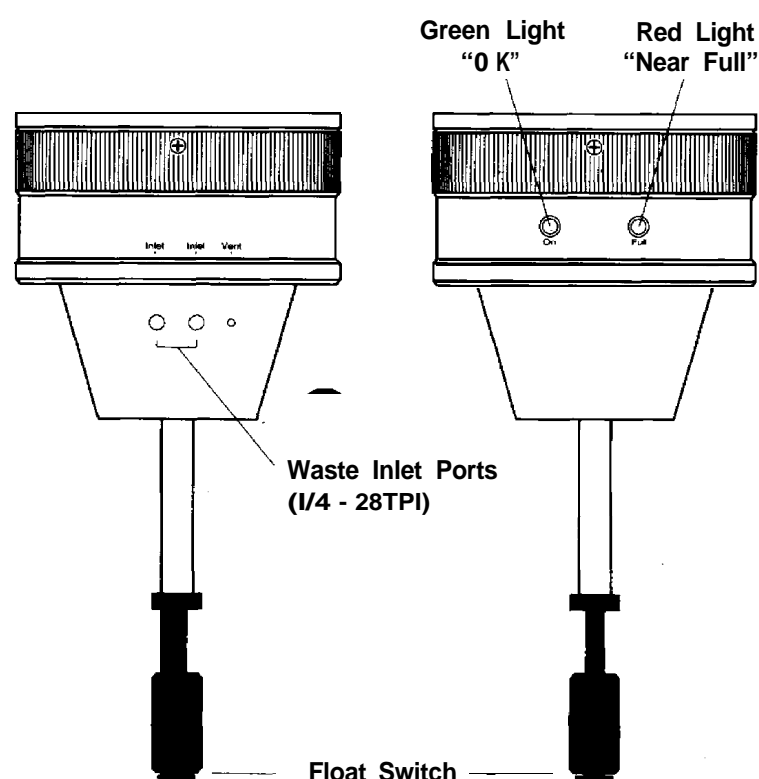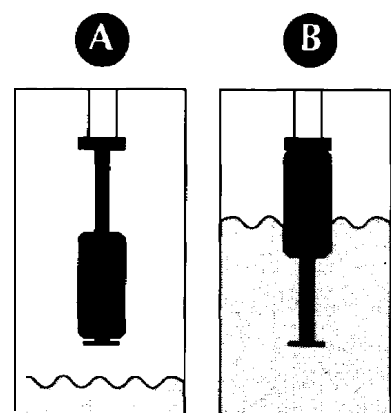

An audible alarm and flashing green and red LED's indicate proper operation and waste bottle condition. Two 1/4-28 TPI ports accept waste solvents via 1/1 6" or tubing.



Level-Minder™ Waste-Bottle Monitor

Part No.	Description
10066	Level-Minder™, complete without bottle
10067	Level-Minder™, complete with safety-coated 2,5 L Amber glass bottle
10068	Level-Minder™, complete with safety-coated 2 L Clear bottle
10069	Replacement Bottle safety-coated 2,5 L Amber bottle
10070	Replacement Bottle safety-coated 2 L Clear bottle
	Larger bottles on request

Distributed by: